
Dry Cleaning of Adhesion Grippers

As an alternative to cleaning adhesion grippers 100015 with water or 70% ethanol, they can also be gently and effectively cleaned using a dry method with a common adhesive tape (e.g. transparent office tape).

Cleaning is recommended when reduced adhesion performance or visible contamination (e.g. particles or dust) is observed on the gripper surface.

SEM image of a microstructured adhesion gripper with thousands of silicone adhesive elements per cm². Adhering contaminants such as particles or dust reduce adhesion and must be removed as required to ensure secure attachment to the bag film.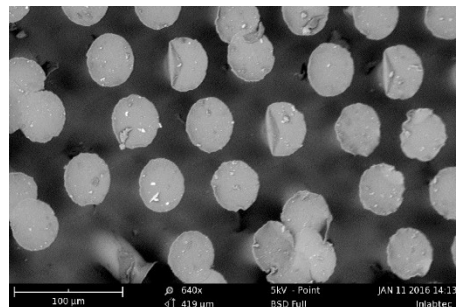

For cleaning and regeneration, the adhesive tape is applied to the surface of the adhesion gripper and lightly pressed down. The tape is then peeled off again. Contaminants adhere to the tape and are removed without the adhesive itself sticking to the silicone material of the adhesion gripper.

The removed contaminants are clearly visible on the peeled-off adhesive tape (image taken with a binocular microscope, 10× magnification).

Notes

- The dry-cleaning procedure may be repeated multiple times if required.
- Use only common transparent adhesive tape.
- Do not use packaging tape, highly adhesive tapes, or aggressive cleaning agents.
- Ensure the gripper surface is dry and free of residues before reuse.

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